


Thin film attachment for ZEM-3

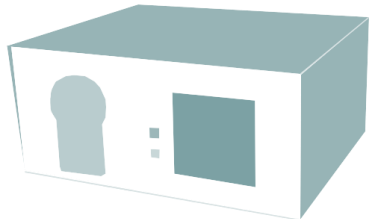

ADVANCE RIKO, Inc.



Alumina screw



Electrode plate



Alumina nut



Washer



Alumina substrate

Procedure 1



(1) Attach electrode plates to both ends of alumina substrate.




Image of (1)



(2) Fix electrode plates to alumina substrate with screw, nut, and washer.



Image of (2) *only one side fixing for stand

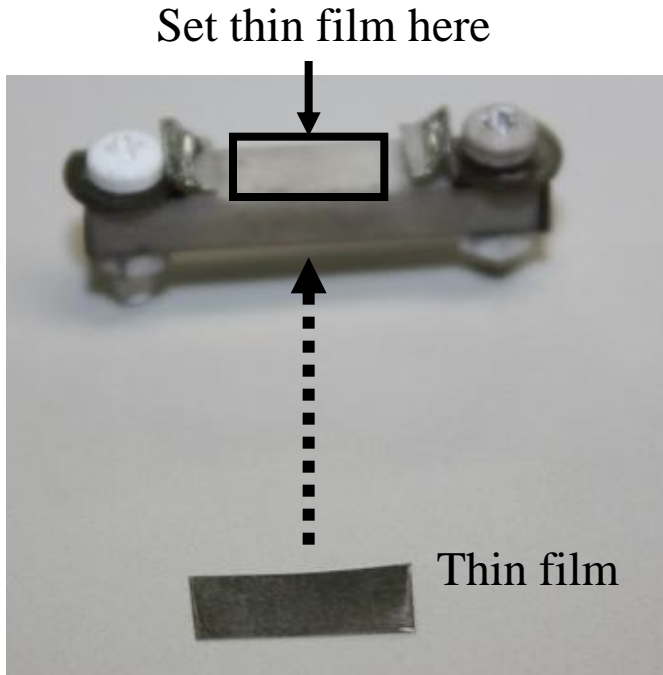

(3) Fix thin film by turning nut.



Image of (3)

* Capable of fixing thin film of various sizes by bending electrode plate.

* If conduction between electrode plate and thin film isn't enough, adhesion should be done with silver paste.

Setting onto ZEM-3

